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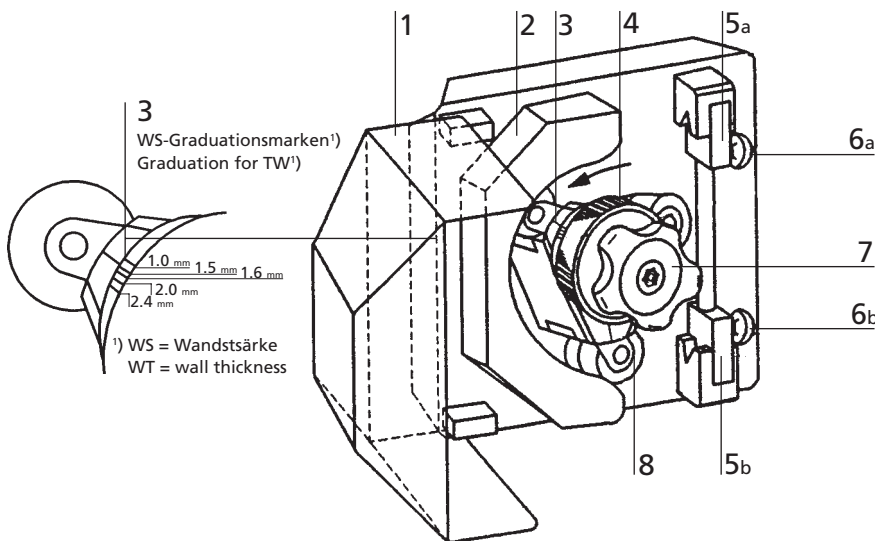
Technical specifications

- Principle: 1-channel tubing pump, with ball bearings
3 stainless steel rollers
adjustable for different tubing types
- Flow-rate: 0.41– 3600 ml/min
with 1.6 mm tubing wall thickness
3.4 – 1500 ml/min
with 2.4 mm tubing wall thickness
- Pressure: max. 2.5 bar (36 psi) with tubing up to
8 mm ID, 1.5 bar (22psi) with larger tubing
- Standard tubing:
WT 1.6 mm ID 1.6 – 11.1 mm
WT 2.4 mm ID 4.8 + 6.4 mm, (WT = wall thickness)

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Adjusting the roller pressure

The knurled ring (4) with the 5 graduation marks allows you to adjust the pressure of the rollers onto the tubing as follows:

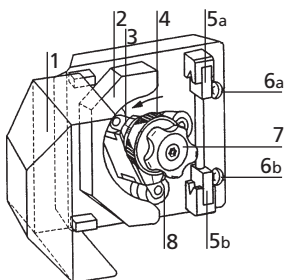
- Open pumphead cover (1)
- Loosen the star grip (7)
→ turn in the counter-clockwise direction
- Turn the knurled ring (4) so that the correct graduation mark (depends on the wall thickness, see figure below) is aligned with the notch in the triangular roller-head (3)
- By simultaneous fastening the star grip (7) you lock the pump-rollers in the position required

The particular advantage of this roller adjustment is that the contact pressure of the 3 rollers occurs concentrically from the roller-head shaft to the contour of the tube-bed. This ensures a regular and constant pressure, independently of the position of the roller-head and the wall thickness of the tubing used. This provides a favorable effect on the pump characteristics and the service-life of the tubing.

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Operating manual for pumphead 380 AD (for the drives MCP oder BVP)

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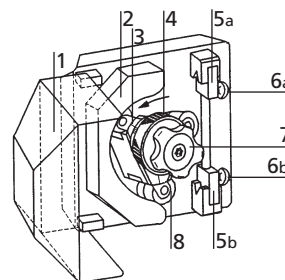
Open the pumphead cover (1)

Load the pump tubing as follows:

- Press the ball handle (6a) down with your thumb
- Insert the tubing with your free hand into the opening of the tubing retainer prism (5a)
- Release the ball handle (6a)
→ the tubing is now clamped in the tubing retainer prism (5a)
- Insert the tubing into the tube-bed (2) by turning the rotor with the pump rollers (8) in the counter-clockwise direction
→ the tubing must be well aligned with the tube-bed (2) and centred by the rollers (8)
- Now lift the ball handle (6b) with your thumb and clamp the tubing in the lower tubing retainer prism (5b)

Start-up procedure

Before switching the drive on, you must close the pumphead cover (1)! Danger of injury!

Before starting the operation of your pump system ensure that the tubing lies correctly in the tube-bed (not twisted) and that the tubing tension is correct. The appropriate tubing tension can be adjusted as follows:

- When pumping in the clockwise direction (*) press down the upper ball handle (6a)
- With your other hand pull slightly the tubing so that a weak tension is achieved
- Release the ball handle (6a)

1) When pumping in the counter-clockwise direction, lift the lower ball handle (6b) and adjust the tubing tension.

→ MCP pumphead identification

ID Code 380A. Must be entered for each of the 4 programs using this pumphead.

Please refer to the drive operating manual.