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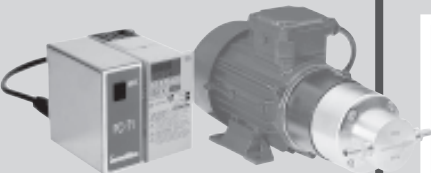
03.10.00 CB/GP

Operating Manual Mode d'emploi

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FC71-MP
ISM 506

Deutsch

Gear pump with variable speed drive

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Contents

Safety precautions	3
Warranty terms	6
Product	8
Operating panel	9
Mains voltage	10
Changing the fuses	11
Before starting the pump	12
Starting the pump	13
Pumping against pressure	14
Before switching the pump off	14
Analog interface (see separate manual for inverter)	14
Pump-head	15
Mounting the pump-head	16
Information on pump-heads	18
Accessories	19
Maintenance, repairs	20
Spare-parts	21
Technical specifications	22

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⚠ Please note

We recommend you to read this operating manual carefully.

We do not take liability for any damage resulting from improper use.

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Safety precautions

Ismatec® gear pumps are designed for pumping applications in laboratories and industry. As such it is assumed that Good Laboratory Practice (GLP) and our following recommendations will be observed.

- The circuit between mains supply and pump has to be earthed.
- The pump must not be operated outside the destined operating and environmental conditions.
- Place the unit in a well ventilated position at least 10 cm away from walls, partitions etc. Ensure that curtains and similar materials do not cover the ventilation slits.
- The pump **must not** be used:
 - for medical applications on human beings
 - in explosion-proof chambers or in the presence of flammable gases or fumes

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⚠ Please note

When operating a gear pump, certain hazards cannot be excluded.

Ismatec SA do not take liability for any damage resulting from the use of an Ismatec® pump.

General risks: Ismatec SA do not admit responsibility for the handling of chemicals.

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Safety precautions

- The pump must be switched off when changing the pump-head.
- The permeability of pump-tubing depends on the material used and pressure conditions. Tubing can also become electro-statically charged. Please be aware of possible hazards when laying tubing in explosion-proof chambers.
- Tubing can tear and burst during operation. If this could cause damage, the necessary safety measures based on the specific situation must be taken.
- Do not manipulate the pump-head before the pump is switched off and disconnected from the mains.
- Before initial start-up, the pump-head must be filled with the pumping liquid or water. Never run the pump-head under dry conditions.

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Warning

Never pump media containing particles.

If necessary, filter the media before starting your application.

Never run the pump-head under dry conditions.

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Safety precautions

- We advise you to never pump liquid containing particles. Even the smallest particles have an abrasive effect, which can ruin the gears in the course of a few seconds.
- Only new fuses, according to the specifications stated on page 11 in this manual, must be used.
- The fuse-holder must not be short-circuited.
- Do not open or remove the housing while the pump is operating.
- Repairs may only be carried out by a skilled person who is aware of the hazard involved.
- For service and repairs carried out by the customer or by third-party companies Ismatec SA denies any responsibility.

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Warranty

From date of delivery: 1 year

In case of any queries, please contact your nearest Ismatec® representative.

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Warranty Terms

We warrant the perfect functioning of the FC71-MP drive, provided it has been installed and operated correctly according to our operating instructions.

If production or material faults can be proved, the defective parts will be repaired or replaced free of charge at our discretion.

The pump-heads are covered by the specific terms of Micropump Ltd. Gears and seals are regarded as consumables and as such they are not covered by the warranty.

The duration of the warranty is not affected by making a claim for warranty service.

Further claims are excluded. Shipping costs for instruments under warranty are charged to the customer.

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Warranty Terms

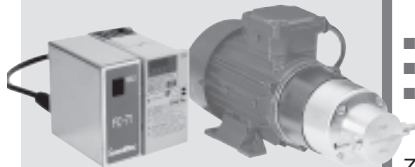
Our warranty becomes invalid in the case of

- improper operation by the user, or if the pump is diverted from its proper use,
- unauthorized modification or misuse by the user or by a third party,
- improper site preparation and maintenance,
- operation outside environmental and electrical specifications for the product,
- use of third-party software, hardware, interfacing or consumables purchased by the user and which do not comply with our specifications,
- damages caused by contamination or leaks due to torn or bursted tubing.

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Product

Package contents

- FC71-MP drive,
incl. magnetic coupling, magnet
and fixing case for the pump-head
- FC71 control unit with frequency
inverter Varispeed 606V7
- Power cord
- Ismatec operating manual
- Instruction manual for the
frequency inverter Varispeed
606V7

Accessories, if ordered:

- Pump-head
- 2 tubing nipples

Please check the package and its contents for transport damage. If you find any signs of damage, please contact your local Ismatec® representative immediately.

Complaints can only be accepted within 8 days from receipt of the goods.

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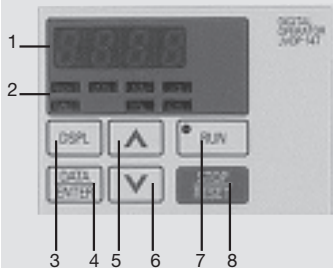
⚠ The drive runs only in a clockwise direction.

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⚠ Please note

For more details see the separate instruction manual for the frequency inverter Varispeed 606V7.

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⚠ Socket/Power cord

Use exclusively the original power cord supplied with the pump. The socket must be earthed (protective conductor contact).

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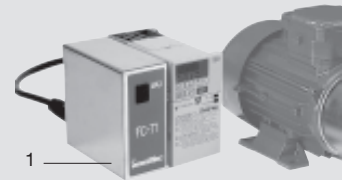
Changing the fuses

- ➔ Switch the pump off, pull out the mains plug.
- 1 Pull out the fuse-holder by opening it with a small screw-driver (size 0).
- 3 Insert 2 new fuses 230 V_{AC}: 2 x 5 A (slow-blow)
 - △ Use always 2 slow-blow fuses of the same type .
- 4 Shut the fuse-holder.

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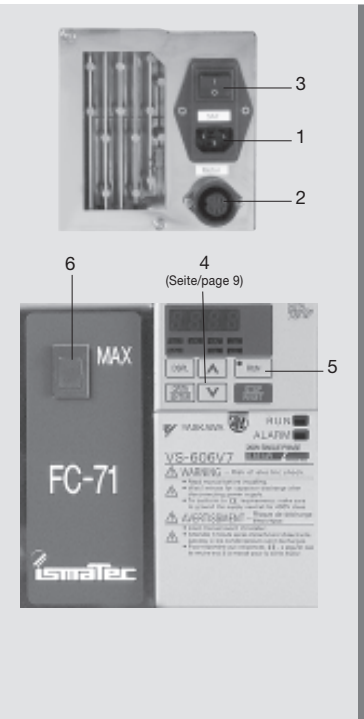
Starting the pump

- 1 Mount the control unit in a vertical position (display and keypad section must be easily accessible) to ensure effective cooling. Mount it on a nonflammable surface (e.g. metal).
- 2 If necessary, mount the pump-head as stated on pages 16–17.
- 3 Before screwing the tubing nipples into the pump-head, we recommend you to wrap PTFE tape around the thread of the nipple. This ensures optimum sealing and helps to protect the thread.
- 4 Mount the tubing and connect it to your system. Fasten the tubing to the nipples by means of the tubing clamps.
- 5 Connect the cable of the motor to the plug on the rear panel of the control unit.

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Starting the pump

⚠ Before starting, fill the pump-head with liquid. The pump-head must never run under dry conditions.

Do not pump media containing particles. If necessary, filter the liquid before use.

- 1 Connect the pump to the mains supply.
- 2 Connect the cable of the motor.
- 3 Press the power supply switch
- 4 Set the required rotation speed (100 = 3450 rpm, max. speed), adjustable in steps of 1%
 ➔ The speed can be adjusted whilst the pump is running.
- 5 Start the pump with the RUN key.
- 6 The MAX- button allows rapid filling (or emptying) of the system whilst the pump is running.

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Pumping against pressure

The max. differential pressure depends on the mounted gear pump-head (see page18).

Before switching the pump off

⚠ If no non-return valve is mounted on the discharge side of the pump, a siphoning may occur at the moment the pump drive is stopped!

If the pump will not be used for some time, it is strongly recommended to thoroughly rinse the pump-head with water.

Analog interface

The frequency inverter features an analog interface.

➔ For further information please refer to the instruction manual of the frequency inverter Varispeed 606V7.

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Pump-head

In general, the FC71-MP gear pump is supplied as a complete unit, consisting of the FC71-MP drive, and if ordered:

- pump-head, incl. magnet
- tubing nipples (2 pcs.)

For the FC71-MP drive, different gear pump-heads with PPS gears are available.

The table on page 18 gives you more detailed information on the pump-heads available.

Before mounting a new pump-head, make sure that the correct magnet is fitted to the drive.

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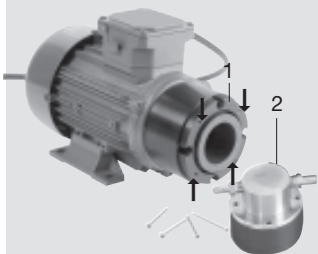
Mounting the pump-head

Fig. 1
For mounting a pump-head, you need a Phillips screwdriver (Size 0).

- 1 Remove all screws from the fixing case of the drive.
- 2 Fasten the pump-head with the 4 long screws that are originally supplied with the drive.

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Mounting the pump-head

Fig. 2

Screw the 2 nipples into the pump-head.

IMPORTANT

Before screwing the tubing nipples into the pump-head, we recommend **you to wrap PTFE tape** around the thread of the nipple. This ensures optimum sealing and helps to protect the thread.

Fig. 3

Screw the pump-head to the fixing-case of the drive. The face of the pump-head with the engraved model number must be directed to the user and the arrow must point to the right.

⚠ Fill the pump-head with liquid before starting to pump. Avoid dry-running of the pump-head.

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Fig. 2

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The casing material is stainless steel (316 ss).
 Range of operating temperature -46°C up to +122°C (-50.8°F up to 251.6°F)
 For temperatures above +54°C (+130°F) consult your Ismatec distributor for proper seal selection. (Please indicate your operation temperature).

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Bestell-Nr. AR 0004

Accessories

Tubing nipple 04 for all gear pump-heads

$\frac{3}{8}$ " thread 18NPT
for tubing i.d. 12.7 mm ($\frac{1}{2}$ "
(always order 2 pcs.)

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Maintenance

After use, we strongly recommend you to thoroughly rinse the pump-head with water.

Provided the FC71-MP drive is operated properly and in compliance with this manual, the moving parts such as gears, etc. are the only parts that are subject to wear and tear.

The Service Kit corresponding to the individual pump-head allows the user to replace the parts that are subject to wear and tear by himself (see page 19).

Repairs

For repairs please send the defective FC71-MP drive to your Ismatec® representative.

Please give information on defect, date of purchase, serial-no., model.

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Spare-parts

After the warranty period your Ismatec® representative will be pleased to send you on request:

- spare-parts
- parts lists
- wiring diagrams

Please give information on defect, date of purchase, serial-no., and model.

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△ Please note

We also recommend you to observe our **Warranty Terms as well as our Terms and Conditions of Sale**.

In case of any queries, please contact your local Ismatec® representative.

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Technical Specifications

Drive

Motor type DC motor
Speed range 0 – 3450 rpm
adjustable in 1% steps

Differential pressure

max. 10.3 bar (149.4 psi)
(depending on pump head
see page 18)

Remote control

via analog interface
(see instruction manual of control
unit)

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Technical Specifications

Operating conditions

Temperature +5 to +40°C,
+41°F to 104°F
Rel. humidity max. 80%
at normal environmental conditions

Dimensions/Weight (Drive)

(drive without pump head)

DxWxH 275x145x180 mm
10.8x5.7x7.1 inches
Weight 8.6 kg / 19 pounds

Commande

DxWxH 160x125x135 mm
6.3x4.9x5.3 inches
Weight 1.9 kg / 4.2 pounds

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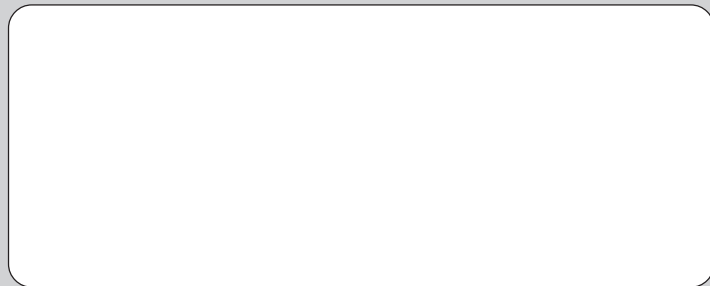
ISMATEC® - Your competent partner for demanding metering and dispensing applications.

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A Unit of IDEX Corporation
Laboratoriumstechnik GmbH