**ISMATEC®**

**IDEX**  
HEALTH & SCIENCE

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03.08.2007 CB/GP

# Operating Manual Mode d'emploi

## Flanschwerkzeug IC 0043

Zubehör (sofern bestellt)

Haltezange IC 0044

Deutsch

## Flanging tool IC 0043

Accessories (if ordered)

Tubing holder IC 0044

English

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Accessoires (si commandés)

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Français

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**Contents**

Safety precautions	4
Warranty terms	6
Product	7
Flanging tool	8
Mains voltage	8
Initial steps of usage	9
Making a flange	10
Accessories	12
Repairs	13
Spare parts	13
Disposal	14
Technical specifications	15

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## Please note

We recommend you to read this operating manual carefully.

When operating an electrically heated device, certain hazards cannot be excluded.

**ISMATEC SA does not take liability for any damage resulting from the use of an ISMATEC® device.**

ISMATEC SA does not admit responsibility for the handling of chemicals.

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## Safety precautions

The ISMATEC® flanging tool has been designed for the use in laboratories. As such it is assumed that Good Laboratory Practice (GLP) and our following recommendations will be observed.

- The circuit between mains supply and flanging tool has to be earthed
- The flanging tool must not be operated outside the destined operating and environmental conditions.
- The flanging tool **must not** be used:
  - in explosion proof chambers or in the presence of flammable gases or fumes
  - in wet zones and areas with splashing water. Note that the housing is ventilated through large slits for thermal reasons. Consequently, the device protection rating is only IP 20.

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## Safety precautions

- Do not touch the heated tip. It gets very hot and there is a risk of getting burned.
- Do not open or remove the housing while the flanging tool is in operation.
- Repairs may only be carried out by a skilled person who is aware of the hazard involved.
- For service and repairs carried out by the customer or by third-party companies ISMATEC SA denies any responsibility.

### **Do not touch the heated tip - risk of burns**

Even several minutes after switching off the flanging tool, the heated tip is still very hot and can cause severe burns.

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## Warranty

### 2 years from date of delivery

We warrant the perfect functioning of our products, provided they have been installed and operated correctly according to our operating instructions.

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## Warranty Terms

If production or material faults can be proved, the defective parts will be repaired or replaced free of charge at our discretion. The duration of the warranty is not affected by making a claim for warranty service. Further claims are excluded. Shipping costs for instruments under warranty are charged to the customer.

### Our warranty becomes invalid in case of

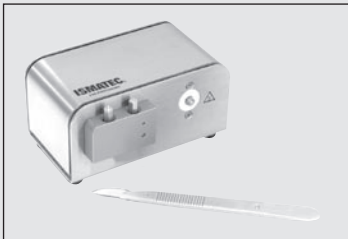
- improper operation by the user, or if the flanging tool is diverted from its proper use,
- unauthorized modification or misuse by the user or by a third party,
- improper site preparation and maintenance,
- operation outside environmental and electrical specifications for the product,
- use of third-party software, hardware, interfacing or consumables purchased by the user and which do not comply with our specifications.

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## Product

### Contents of the package

- Flanging tool IC 0043
- Power cord
- 3 different heated tips:  
Small for tubing i.d. 0.3 mm  
Medium for tubing i.d. 0.5/0.8 mm  
Large for tubing i.d. 1.5/2.0 mm
- 1 scalpel
- Operating manual

Accessories (if ordered)  
Tubing holder IC 0044

Please check the package and its contents for transport damage. If you find any signs of damage, please contact your local ISMATEC® representative immediately.

Complaints can only be accepted within 8 days from receipt of the goods.

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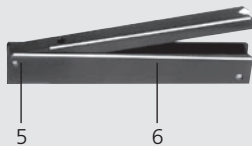
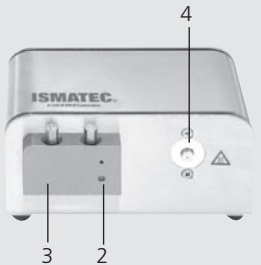
## Flanging tool

- 1 Mains socket
- 2 Hole for adjusting the correct length of the tubing tip
- 3 Plate
- 4 Heated tip
- 5 Opening for inserting the PTFE tubing
- 6 Tubing holder

## ⚡ Mains voltage

100 – 240 V<sub>AC</sub> 47 – 63 Hz

### ⚠ Socket/Power cord

Use exclusively the original power cord supplied with the flanging tool. The socket must be earthed (protective conductor contact).

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## Initial steps of usage

- 1 Select the heated tip and screw it into the flanging device. (choice of the tip: s. p. 7)  
The rear end of the tubing holder serves as a wrench.
- 2 Connect the flanging tool to the mains using the original power cord.
- 3 Wait about 10 to 15 minutes until the device has reached the operating temperature.

### Do not touch the heated tip - risk of burns

Even several minutes after switching off the flanging tool, the heated tip is still very hot and can cause severe burns.

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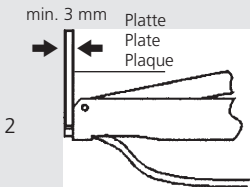
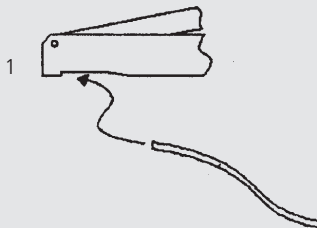
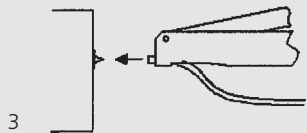
## Making a flange


- 1 When cutting the PTFE tubing ensure that the tubing ends are cut precisely at right angles. Otherwise the flange surface will not be tight.
  - ➔ Consider that after producing a flange, the length of the tubing will be shortened by 2 mm.
  - ➔ Use a scalpel in order to get a precise cut.
- 2 Push the PTFE tubing through the hole in the tubing holder and insert it in the corresponding groove:
  - Tubing 0.3 / 0.5 / 0.8 i.d. mm: small groove
  - Tubing 1.5 / 2.0 i.d. mm: large groove.
  - ➔ The tubing must jut out by at least 3 mm. Use the corresponding hole on the plate of the flanging tool in order to determine the correct length of the protruding tubing:
    - Tubing 0.5 / 0.8 mm i.d.: small hole
    - Tubing 0.3 / 1.5 / 2.0 mm i.d.: large hole.

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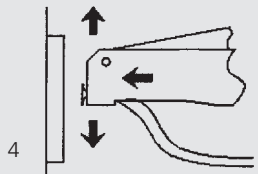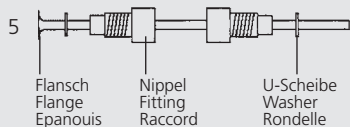
## Making a flange

- 3 Compress the tubing holder and press the tubing at right angles onto the heated tip. At the same time you slowly rotate repeatedly the tubing to the left and the right.
- 4 Press the flange at right angles against the plate and move it up and down about 4 to 5 times, until the flange has cooled down and the material is hard again.
- 5 Before starting to produce the second flange, both fittings and washers must be pushed onto the tubing, as illustrated on the left. For the second flange repeat paragraphs 1 to 4.

6 Completed tubing connection.

**Do not touch the heated tip! Risk of getting burned!**

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Tubing i.d. 0.5 and 0.8 mm IS 1076

Tubing i.d. 1.5 mm IS 1075

→ For fittings, tubing, etc.,  
please refer to our catalog.

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## Repairs

For repairs we ask you to send the complete flanging tool with detailed description of the failure to your ISMATEC® distributor. Please use the original ISMATEC® packing or one of equal quality.

## Replacement parts

For repairs you intend to carry out by yourself (out of the warranty period), you can ask from your ISMATEC® distributor:

- Parts lists
- Replacement parts
- Wiring diagrams

Please give information on defect, date of purchase, serial No., and model.

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## Disposal

Please retain packing materials until the product warranty ends. Afterwards please discard packing materials in an environment-friendly manner according to local regulations.

Once the useful life of the product has ended, please ensure proper disposal according to local laws. Plastic and electronic components should be disposed of at a recycling facility. Please refer to local regulations regarding proper disposal.

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## Please note

We also recommend you to observe our Warranty Terms as well as our Terms and Conditions of Sale.

In case of any queries, please contact your local ISMATEC® representative.

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Temperature +41°F to 104°F

Rel. humidity max. 80 %

– not condensing, at normal laboratory conditions

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Ask for the latest  
ISMATEC catalogue.

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ISMATEC® - Your competent partner for demanding metering and dispensing applications.

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Laboratoriumstechnik GmbH