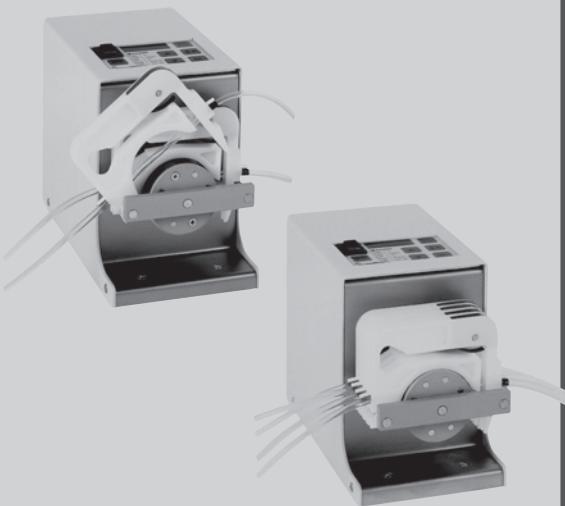
REGLO *Digital*

Schlauch- pumpe

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Schutzgrad IP 30

2 Kanäle

ISM 831	6 Rollen
ISM 832	8 Rollen
ISM 596	12 Rollen

4 Kanäle

ISM 833	6 Rollen
ISM 834	8 Rollen
ISM 597	12 Rollen

Tubing pump

Microprocessor
controlled

Protection rating IP 30

2 channels

ISM 831	6 rollers
ISM 832	8 rollers
ISM 596	12 rollers

4 channels

ISM 833	6 rollers
ISM 834	8 rollers
ISM 597	12 rollers

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Classe de protection IP 30

2 canaux

ISM 831	6 galets
ISM 832	8 galets
ISM 596	12 galets

4 canaux

ISM 833	6 galets
ISM 834	8 galets
ISM 597	12 galets

Deutsch

English

Français

**ISMATEC**[®]

IDEX
IDEX CORPORATION

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⚠ Please note

We recommend you to read this operating manual carefully.

When operating a pump, certain hazards cannot be excluded.

ISMATEC SA does not take liability for any damage resulting from the use of an ISMATEC® pump.

ISMATEC SA does not admit responsibility for the handling of chemicals.

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Safety precautions

ISMATEC® Pumps are designed for pumping applications in laboratories and industry. As such it is assumed that Good Laboratory Practice (GLP) and our following recommendations will be observed.

- The circuit between mains supply and pump has to be earthed.
- The pump must not be operated outside the destined operating and environmental conditions.
- The pump must not be used:
 - for medical applications on human beings
 - in explosion proof chambers or in the presence of flammable gases or fumes
- The pump must be switched off when pump-heads, cassettes or tubing are inserted or changed.
- The permeability of tubing depends on the material used and pressure conditions. Tubing can also become electro-statically charged. Please be aware of possible hazards when laying tubing in explosion-proof chambers.
- Pump-heads consist of revolving parts. Therefore, the pump must not be operated before the cassettes are fully snapped-in.

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Safety precautions

- Do not manipulate the pump-head before the pump is switched off and disconnected from the mains supply.
- Be particularly cautious that no parts of your body such as fingers, long hair, etc. or jewellery, or loose objects such as cables or tubing, etc. can be trapped by the revolving rotor.
- Tubing can tear and burst during operation. If this could cause damage, the necessary safety measures based on the specific situation must be taken. When pumping corrosive media we recommend to place the pump in a collecting basin.
- Only new fuses, according to the specifications stated on page 9 in this manual, must be used.
- The fuse-holder must not be short-circuited.
- Do not open or remove the housing while the pump is operating.
- Repairs may only be carried out by a skilled person who is aware of the hazard involved.
- For service and repairs carried out by the customer or by third-party companies ISMATEC SA denies any responsibility.

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Warranty

For all parts manufactured by
ISMATEC®
All other parts,
excluding consumables: **1 year**

In case of any queries, please
contact your ISMATEC®
representative.

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Warranty terms

We warrant the perfect
functioning of our products,
provided they have been installed
and operated correctly according
to our operating instructions.
If production or material faults
can be proved, the defective
parts will be repaired or replaced
free of charge at our discretion.

A pump must be returned in the
original ISMATEC® packing or in a
packet of equal quality. The
duration of the warranty is not
affected by making a claim for
warranty service.

Further claims are excluded.

Shipping costs are charged to the
customer.

Our warranty becomes invalid in the case of:

- improper operation by the
user, or if the pump is diverted
from its proper use
- unauthorized modification or
misuse by the user or by a
third party
- improper site preparation and
maintenance
- operation outside of the
environmental and electrical
specifications for the product
- use of third-party software,
hardware, accessories or
consumables purchased by the
user and which do not comply
with our specifications.

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Contents of the package

- Pump REGLO *Digital* (type as ordered)
- 2 or 4 cassettes
- 1 power cord with country specific mains plug
- Operating manual

Please check the package

and its contents for transport damage. If you find any signs of damage, please contact your local ISMATEC® representative immediately.

Complaints can only be accepted within 8 days from receipt of the goods.

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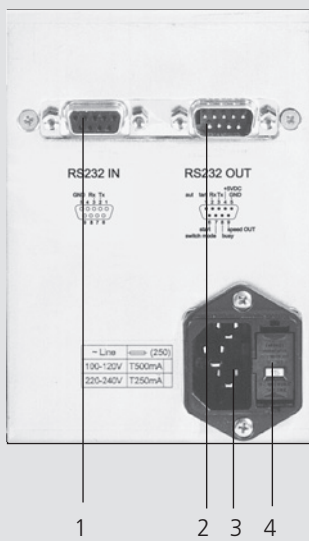
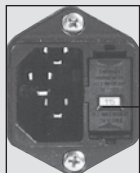
Rear panel

- 1 RS232 IN (female) for operation controlling via PC
- 2 RS232 OUT (male) for cascading up to 8 pumps

Digital input (TTL level)
Run/Stop, autostart

Analog output

speed output:

- 0–8 kHz (2-channel version)
- 0–5 kHz (4-channel version)

- 3 Mains socket
- 4 Fuse-holder with voltage selector 115/230V

⚡ Mains voltage

Mains voltage	Fuse rating
220-240 V _{AC}	2 x 250 mA *)
110-120 V _{AC}	2 x 500 mA *)

*) slow-blow type

⚠ Socket/Power cord

Use exclusively the originally supplied power cord.
The socket must be earthed (protective conductor contact).

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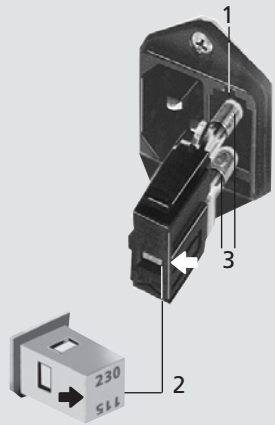
Voltage selector and
window in the fuse-holder

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Voltage setting 115V / 230V and changing the fuses

➔ Switch the pump off,
pull out the mains plug.

- 1 Pull out the fuse-holder by opening it with a small screw-driver (size 0).
- 2 Take out the voltage selector plate. Turn it and re-insert it into the fuse-holder so that the required voltage rating is facing the window of the fuse-holder.
- 3 Insert 2 new fuses
230 V_{AC}: 2 x 250 mA (slow-blow)
115 V_{AC}: 2 x 500 mA (slow-blow)

⚠ Use always 2 slow-blow fuses of the same type complying with the local mains voltage.
- 4 Shut the fuse-holder. The voltage rating is visible in the window.

⚠ Before starting-up

Check if the voltage setting visible in the window of the fuse-holder complies with your local mains voltage.

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Operating panel

- 1 On/Off switch
- 2 Digital LED display
- 3 Run/Stop
- 4 Double function:
 - Rotation direction
 - Reset
- 5 Calibration / Saving a set value
- 6 Reduce value
- 7 Increase value
- 8 MODE key
 - for selecting operating mode
 - Escape key for value setting

9 LEDs for active operating mode

- **PUMP • Flow rate**
Pumping by flow rate (ml/min) (also by speed if rpm is chosen for this mode).
➔ In 'Tube • mm I.D.' you can switch to rpm (see page 14).
- **DISP • Volume**
Dispensing by volume (ml) (or by roller steps – if rpm is chosen for this mode – see page 14).
- **DISP • Time**
Dispensing by time
- **PAUSE • Time**
Pause time for dispensing at intervals
- **TUBE • mm I.D.**
Tubing i.d. selection and changing to rpm

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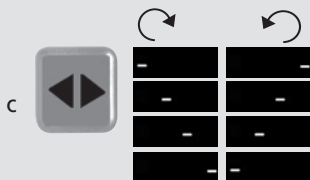
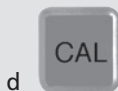
Control keys

a RUN/STOP

- ➔ Starts and stops the pump
- ➔ Starts, interrupts and continues a dispensing cycle

b MODE

- Changes between operating modes (see p. 10)
- Escape key for value setting

c Rotation direction

Changes the rotation direction.

By changing the rotation direction or starting the pump, the rotation direction is indicated in the display with a running hyphen.

Reset key for:

- resetting an interrupted dispensing cycle
- resetting a titration volume to ZERO
- Default key for value resetting, e.g. for resetting a flow rate or volume calibration to its default value (see p. 20)

d CAL

Multi-function key for:

- calibrating the flow rate and dispensing volume
- max. speed (when pump is running)
- saving a set value

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Start-up information

After switching on the power supply switch, the following values are displayed:

- 1 LED test »8.8.8.8.«
- 2 Firmware version
- 3 Pump type (number of channels, number of rollers)
- 4 Roller back-steps
➔ are only displayed if at least 1 or more roller back-steps are entered
- 5 Displays the operating mode last used

Resetting the parameters to the default values

- 1 Power switch »OFF«
- 2 Push the **MODE key** whilst switching the pump on

The following parameters are reset to the following default values:

Mode: PUMP • Flow rate
Rotation direction: clockwise

Flow rate:*
depends on pre-set tubing i.d.

Dispensing volume:*
depends on pre-set tubing i.d.

Dispensing time:
4.5 sec

Pause time:
2.0 sec

***is based on 40.0 rpm**

Dispensing cycles: 0 (infinite)
Roller back-steps: 0
Calibration: Default calibration

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Ersatz-Kassetten

MS/CA Click'n'go	
Bestell-Nr.	IS 3510
MS/CA Anpresshebel*	
Bestell-Nr.	IS 0649

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Inserting the tubing

△ First enter the tube i.d. as described on page 14.

- ➔ Switch the pump off
- Remove the cassette by slightly pressing the fixing-tongue and lifting it simultaneously
- Insert the 3-stop color-coded tubing with one stopper into the cassette
- Let the tubing hang down (prevent it from being twisted)
- Insert the tubing with the second stopper at the other end of the cassette
- Reinsert the cassette into the roller-head

- ➔ Switch the pump on

➔ When the pump is idle, release all cassettes at the fixing-tongue (2).

➔ Flow rate chart see page 37

Spare-cassettes

MS/CA Click'n'go	
Order-No.	IS 3510
MS/CA pressure lever*	
Order No.	IS 0649

* The cassettes with pressure lever are available on request. This type of cassette may provide better results both for tubing Tygon MH and at elevated differential pressure conditions.

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


Cassettes de rechange

MS/CA Click'n'go	
No de commande	IS 3510
MS/CA levier de pression*	
No de commande	IS 0649

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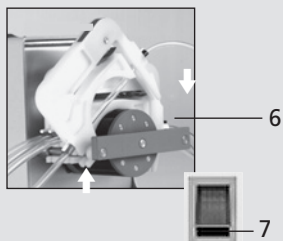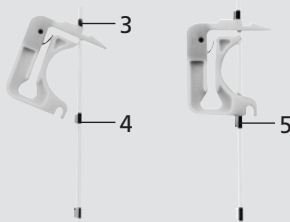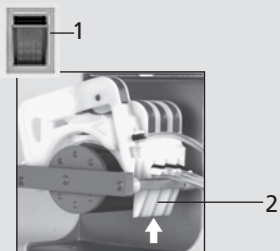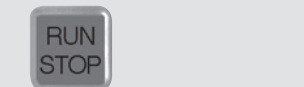
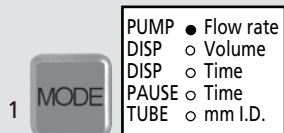
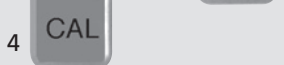
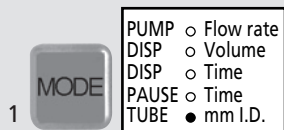
3-stop color-coded tubing

Please refer to our detailed tubing documentation.

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Setting the tube size (i.d.)

- 1 Change MODE to TUBE • mm I.D.
- 2 Press CAL-key (display is blinking)
- 3 Enter the required tube i.d. with ▲▼ keys
- 4 Confirm value with CAL-key
 ➔ Change to the required operating mode by using the MODE key

⚠ In mode TUBE • mm I.D. you can also select 'rpm' (speed), which has the following consequences:

- In mode PUMP • Flow rate the revs per minute (rpm) are indicated.
- In mode DISP • Volume you can read the number of roller steps instead of the volume (ml).

Pumping by flow rate

- 1 Change mode to PUMP • Flow rate
- 2 Enter the required flow rate with ▲▼ keys (displayed in ml/min) Start pump with RUN/STOP
 ➔ The flow rate can also be adjusted while the pump is running.
 ➔ For an accurate flow rate we recommend to calibrate the pump (see page 15).
- 3 For fast filling or emptying the system, maintain pressure on the CAL key (while the pump is running).

⚠ If you select 'rpm' in the mode Tube • mm I.D., the value displayed in PUMP • Flow rate will be the speed (rpm) instead of the flow rate (see paragraph above).

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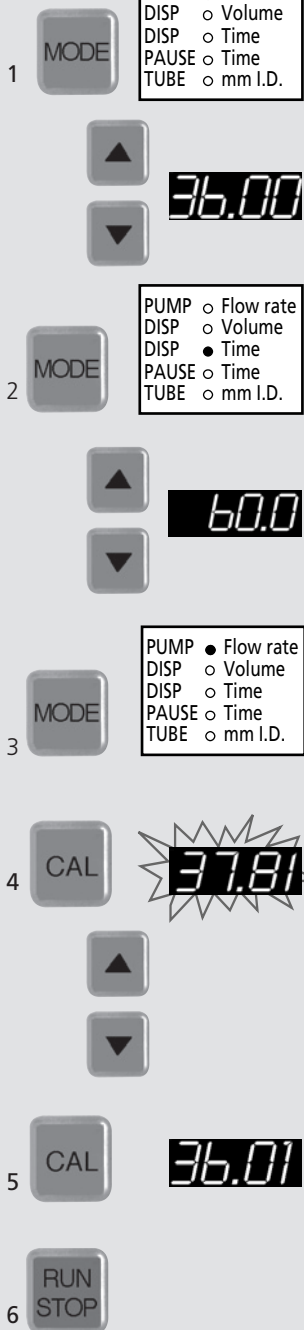
Calibrating the flow rate

- Change mode to PUMP • Flow rate
Enter the required flow rate with the ▲▼ keys
➔ Check whether the correct tubing i.d. is selected in the mode Tube • I.D. mm?
- Change mode to DISP • Time
– Enter 60 seconds by using the ▲▼ keys
– Start pump with RUN/STOP
➔ Pump stops automatically after 60 seconds
➔ Measure the dispensed liquid by volume or weight.

Calibrate the ascertained value as follows:

- Return to mode PUMP • Flow rate
- Press the CAL-button (displayed value blinks)
Enter the measured value (balance) with the ▲▼ keys
- Confirm with the CAL-key (the flow rate setting is readjusted to the initially entered set point as close as possible).
The rotation speed has a resolution of $\frac{1}{10} \text{ rpm}$.
After calibration, very often the initially set flow rate can only be reached approximately.

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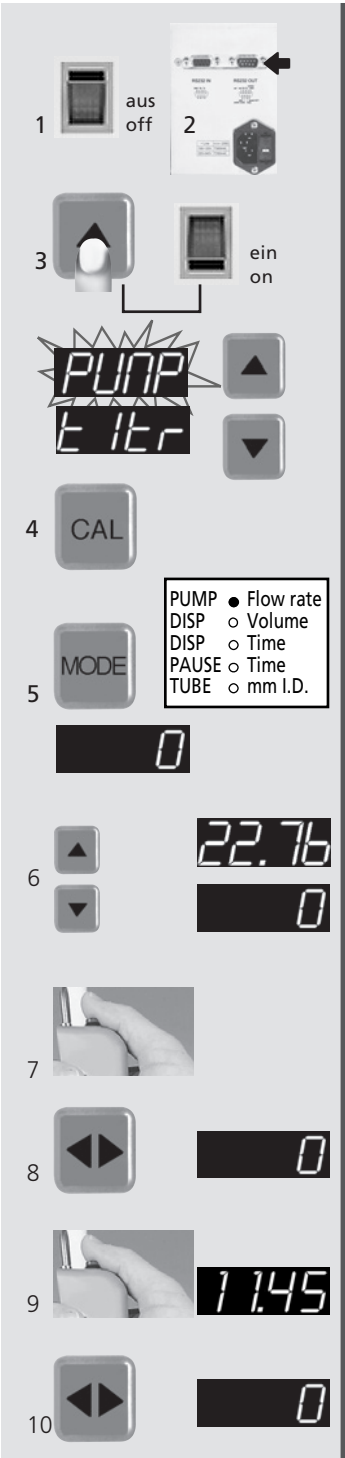
Titration

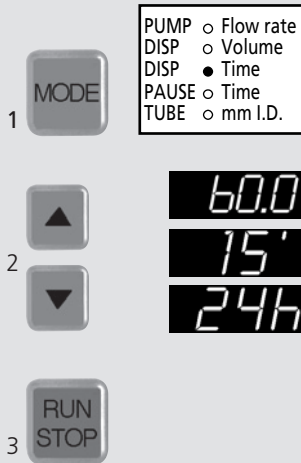
➔ For an accurate titration we recommend you to calibrate the pump(see page 19).

- 1 Switch the pump off.
- 2 Connect the titration handle to RS232 OUT.
- 3 Keep ▲ key pressed and switch power 'ON' (displayed value blinks)
 - ➔ Change to 'titr' with the ▲▼ key
- 4 Confirm with the CAL-key
- 5 Change mode to PUMP • Flow rate
 - ➔ '0' appears in the display (titration volume).
- 6 Set the required flow rate with the ▲▼ keys. (is indicated in ml/min)
 - ➔ Afterwards the reading in the display returns to '0'.
- 7 Start the pump with the titration handle and fill the tubing line.
 - ➔ Maintain pressure on the push button.
- 8 Reset the titration volume to '0' by using the ◀ key.
- 9 Start the pump with the titration handle and carry out the titration.
 - ➔ The push button must be pressed down during the titration.
 - ➔ The titrated volume is shown in the display
- 10 The volume can be reset to '0' by using the ▶ key.

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Please note

- A dispensing cycle can be interrupted by pressing the RUN/STOP key.
- the remaining dispensing time blinks
 - by pressing the RUN/STOP key again, the dispensing cycle is continued
 - by pressing the **◆**-key an interrupted dispensing cycle can be stopped for good

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Dispensing by time

The dispensing time can be entered from 0.1s to 999h.

- 1 Change mode to DISP • Time
- 2 Enter the required dispensing time with the **▲▼** keys
 ➔ The display accelerates when pressure on the **▲** or **▼** key is maintained.
 Seconds: 0.1–899.9 (in 0.1s steps)
 Minutes: 15'–899' (in 1min steps)
 Hours: 15h–999h (in 1h steps)

3 Start pump with RUN/STOP

➔ With the **▲▼** keys the dispensing time can be changed while the dispensing cycle is in process.

➔ During the dispensing process the flow rate (PUMP • Flow rate) can be retrieved by pressing the MODE-key. The flow rate can be changed during the dispensing process by using the **▲▼** keys.

➔ Very short dispensing times can result in non-repeatable dispensing volumes.

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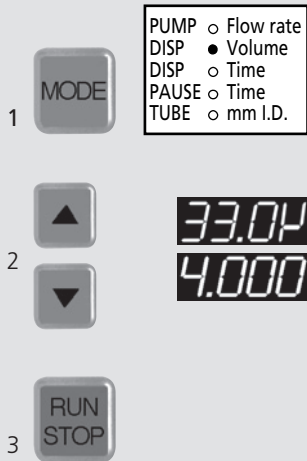
Please note

A dispensing cycle can be interrupted by pressing the RUN/STOP key.

- the volume dispensed up to now is blinking
- by pressing the RUN/STOP key again, the dispensing cycle is continued
- by pressing the key the dispensing cycle is stopped for good

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Dispensing by volume

- 1 Change mode to DISP • Volume
- 2 Use the keys for entering the required dispensing volume. Two display readings are available:
Microlitre: e.g. 12.5 µl = **12.5µ**
Millilitre: e.g. 230.5 ml = **230.5**
➔ For dispensing accurately we recommend you to calibrate the pump (see p. 19).
- 3 Start pump with RUN/STOP. The pump reduces the rotation speed shortly before the end of the dispensing cycle providing controllable and drop-precise dispensing cycles.

➔ The dispensing speed can be adjusted in the mode PUMP • Flow rate.

➔ With the keys the dispensing volume can also be changed even during the dispensing process. The newly entered set point appears shortly in the display.

➔ During the dispensing process the flow rate (PUMP • Flow rate) can be retrieved by pressing the MODE-key.

➔ The flow rate can be changed even during the dispensing process by using the keys .

If you select 'rpm' in the mode Tube • mm I.D. , the number of roller steps will be indicated in DISP • Volume (see page 14)

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Calibrating the volume

- 1 Change MODE to DISP • Volume
➔ Use the ▲▼ keys for entering the required dispensing volume
- 2 Start with RUN/STOP
➔ The pump reduces the rotation speed shortly before the end of the dispensing cycle providing controllable and drop-precise dispensing volumes
➔ Measure the dispensed liquid by volume or weight
- 3 Stay in the same mode (DISP • Volume) and press the CAL-key (display blinks)
- 4 Enter the effectively dispensed volume with the ▲▼ keys
- 5 Save setting with CAL-key
➔ The calibrated volume is adjusted to the initially entered set point as close as possible. (The volume per roller step determines the smallest calibration unit. A dispensing volume is always the multiple of this unit. Very often, the initially set volume can only be reached approximately).
- 6 Start pump with RUN/STOP

➔ Depending on the application conditions repeated calibration may be necessary.
➔ The smallest possible dispensing volume corresponds to the volume of one roller step.

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Default calibration of flow rate

- 1 Change mode to PUMP • Flow rate.
 - 2 Press the CAL button (displayed value blinks)
 - 3 Press the ◀▶ key. (Default value blinks).
 - 4 Confirm with the CAL button.
- ➔ At the same time, the volume is also reset to the default value.

Default calibration of volume

- 1 Change mode to DISP • Volume
 - 2 Press the CAL button (displayed value blinks)
 - 3 Press the ◀▶ key. (Default value blinks).
 - 4 Confirm with the CAL button
- ➔ At the same time, the flow rate is also reset to the default value.

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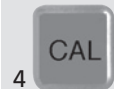
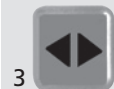
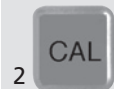
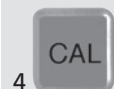
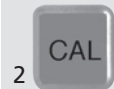
Dispensing by volume within a pre-set time

1 Change to mode DISP • Volume . Enter the required volume (in µl or ml) with the ▲▼ keys.

2 Change to mode DISP • Time. Enter the required time with the ▲▼ keys (0.1s – 999h) (see page 17)

3 Push the MODE-button once
➔ Both LEDs light up
➔ 1111 and 9999 indicate an error message (see below)

4 Start pump with RUN/STOP
➔ The pump dispenses the required volume within the pre-set time.

a Changes / Corrections
For changing the dispensing time or dispensing volume press the ◀▶ key.
➔ The pump returns directly to MODE DISP • Volume
➔ Repeat steps 1 to 3

b Error message
LEDs are blinking alternately
9999 = volume too large
time too short
1111 = volume too small
time too long

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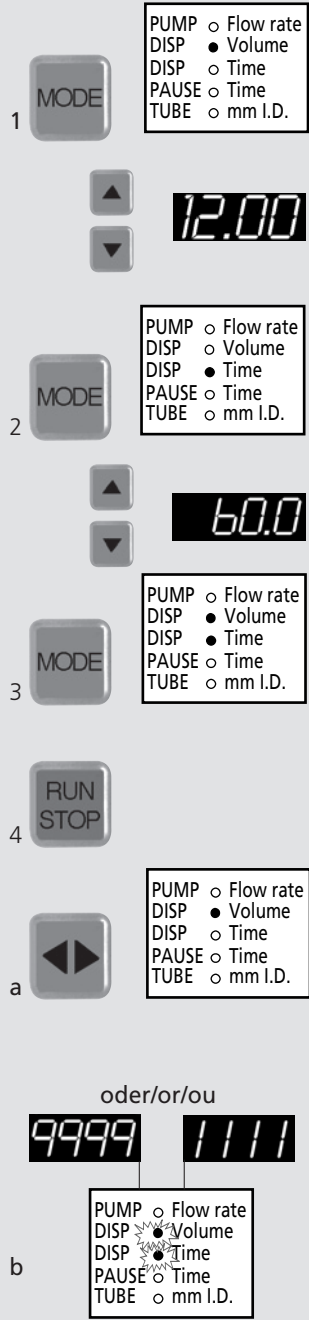
Intermittent dispensing (by time)

Intermittent dispensing by time with a pre-set pause time

- 1 Go to mode PAUSE • Time
Enter the pause time (between 0.1s and 999h) with ▲▼ keys.
- 2 Press MODE key until both LEDs for PAUSE and DISP • Time light up simultaneously.
➔ Enter the dispensing time (0.1s to 999h) with ▲▼ keys
- 3 Start with the RUN/STOP key.
➔ The pump dispenses now at intervals.
➔ Active mode blinks.
- 4 Pushing the RUN/STOP key again interrupts the dispensing process or pause period.
➔ The display blinks.

The user can now proceed as follows:

- ➔ By pressing the RUN/STOP key again, the dispensing or pause period is continued.
- ➔ By pressing the  key the pump is stopped for good.

- ➔ With the ▲▼ keys the dispensing or pause time can be changed even during the dispensing process. The newly entered set point appears shortly on the display.
➔ During the dispensing process the flow rate can be retrieved with the MODE button and adjusted with the ▲▼ keys.

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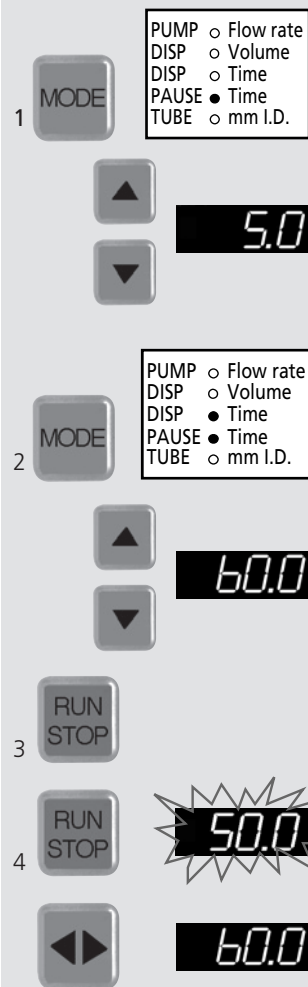
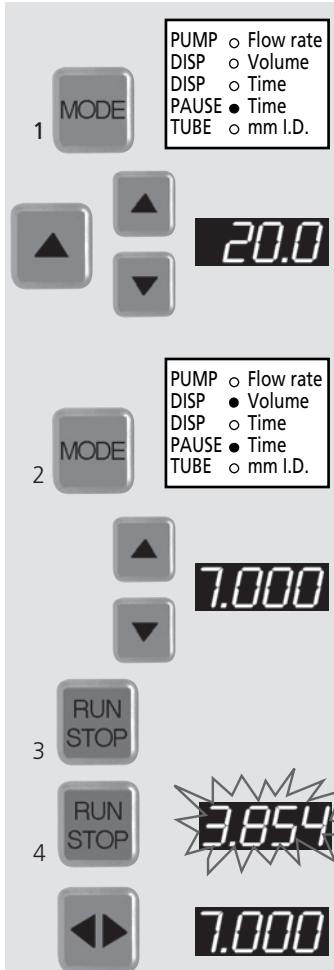
Please note

For entering the number of dispensing cycles see page 24.

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PUMP ○ Flow rate
DISP ○ Volume
DISP ○ Time
PAUSE ● Time
TUBE ○ mm I.D.

PUMP ○ Flow rate
DISP ● Volume
DISP ○ Time
PAUSE ● Time
TUBE ○ mm I.D.

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Intermittent dispensing (by volume)

Intermittent dispensing by volume with a pre-set pause time

- Go to mode PAUSE • Time. Enter the pause time (between 0.1s and 999h) with ▲▼ keys.
 - Press MODE key until both LEDs for PAUSE and DISP • Volume light up simultaneously. ➔ Enter the dispensing volume with the ▲▼ keys.
 - Start with the RUN/STOP key. ➔ The pump dispenses now at intervals. ➔ Active mode blinks.
 - Pushing the RUN/STOP key again interrupts the dispensing process or pause period. The user can now proceed as follows:
 - By pressing the RUN/STOP key again, the dispensing or pause period is continued.
 - By pressing the ◀ key the pump is stopped for good.
- ➔ With the ▲▼ keys the dispensing volume or pause time can be changed even during the dispensing process. The newly entered set point appears shortly on the display. ➔ During the dispensing process the flow rate can be retrieved with the MODE button and adjusted with the ▲▼ keys .

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Number of dispensing cycles

For intermittent dispensing (by time or volume), the number of dispensing cycles can be entered.

- 1 Power switch »OFF«
- 2 Push the ▼ key whilst switching the pump on
- 3 Enter the number of dispensing cycles with the ▲▼ keys.
➔ 0 = continuous dispensing
- 4 Confirm with the CAL-key
- 5 Change into one of the two modes for dispensing at intervals
– DISP•Time / PAUSE•Time or
– DISP•Volume / PAUSE•Time
- 6 Start the pump with the RUN/STOP-key

➔ During the pauses the display shows the remaining number of dispensing cycles.

➔ During long pauses, the remaining interval time can be checked with the CAL-key.

➔ At any time, the dispensing process or pause period can be interrupted with the RUN/STOP key.

➔ By pressing the RUN/STOP key again the dispensing process or pause is resumed with the remaining amount of dispensing cycles.

➔ An interrupted dispensing cycle can be stopped for good by using the ◀ key.

➔ The number of dispensing cycles remains saved even when the pump is switched off.

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Drip-free dispensing

With programmable roller back-steps (1 to 100 steps)

- 1 Switch the pump off
- 2 Maintain pressure on key and switch the pump on again
- 3 Display blinks
Enter the number of roller back-steps with the keys
- 4 Save input with the CAL key

Roller back-steps are active in the following modes:

- a Dispensing by volume (see page 18)
- b Dispensing a volume within a pre-set time (see page 21)
- c Intermittent dispensing by volume (see page 23)

Reject the first dispensing step. Afterwards, the roller back-steps are compensated.

Deleting roller back-steps

- Repeat step 1 and 2
- Enter figure 0 with the keys
- Save input by pressing the CAL key

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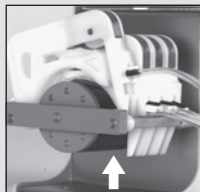
Pumping against pressure

The REGLO *Digital* can be used for continuous duty at a differential pressure of max. 1.0 bar (smaller tube sizes and/or lever type cassettes allow higher pressures)

When the pump is idle

When the pump is idle, we recommend you to release the tubing from pressure. Releasing the cassette on the right side is sufficient. This helps to protect the tubing from unnecessary strain and prolongs its service-life.

⚠ Syphoning effect

When the tubing is released from squeezing, the fluid can flow back to the reservoir.

Overcurrent protector

The REGLO *Digital* drive features an overload protector. When an overload condition occurs, it is indicated in the display by the letters 'OL' and the pump is stopped.

Whenever this situation occurs, **the pump must be switched off immediately**. Before the pump is re-started, it is most important to check the reason for the overload (eg. too high differential pressure).

➔ Only when the cause of the overload has been detected and the failure corrected accordingly may the pump be started again.

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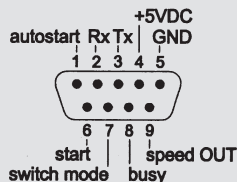
⚠ Please note

The analog interface is integrated in the port 'RS232 OUT' and can be used for remote control (e.g. by foot-switch).

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Analog interface

Pin 1: autostart

If this input is connected to the ground, the pump will automatically start in the operating mode set before the pump was switched off.

Pin 4: +5VDC

For connecting an external control unit (current <100mA).

Pin 5: GND

Point of reference for all signals described.

Pin 6: start

Connection for an external trigger device (hand dispenser, foot-switch, etc.) For operating modes, refer to pin 7.

Pin 7: switch mode

- Input open:
Pump starts or stops when the external trigger device is applied
- Connection to ground:
Pump is running as long as the external trigger device is pressed

Pin 8: busy

Informs about pump status.

Pumpe in operation

= output +5V

Pumpe in stand-by

= output on ground

(The output resistance is 4.7 kΩ)

Pin 9: speed OUT

A square-wave signal with a frequency proportional to the speed (rpm) is at your disposal: 3000 pulses are generated per one rotor revolution.

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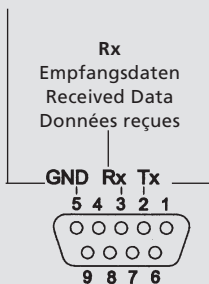
⚠ Please note

When switching on the REGLO *Digital* pump, the serial interface will only respond after a delay of 3 seconds.

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Serial interface

RS232 IN (female)

A 9-pin D-socket is available on the rear panel of the pump.

Setting the pump address

Each command string begins with the pump address (1–8). It is factory-set at 1 and can be changed with @. The address will be stored permanently (even after the pump has been switched off). The assignment of the address enables the user to control up to 8 ISMATEC® drives via one interface.

Structure of the commands

The address is followed by a character. Some commands have an additional parameter which always consists of 4 or 5 figures.

The command string is **completed by** a carriage return (ASCII 13). The pump confirms most of the commands with an asterisk *. Yes/No inquiries are answered by + (yes) or – (no).

Multi-digit replies are concluded by (cr) ASCII 13 and (lf) ASCII 10.

Incorrect command strings are answered by #.

Numerical values are confirmed as 3 to 5-digit figures. Four of the five digits are numerals, one digit is either a decimal point or a preceding blank space.

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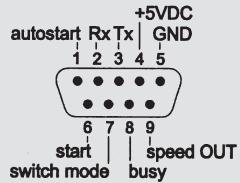
RS232 OUT (male)

Pin 2: RS232 Rx

Is used for connecting additional pumps together with pin 3/5.

Pin 3: RS232 Tx

Is used for connecting additional pumps together with pin 2/5.

Pin 5: GND

Point of reference for all signals described.

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** Roller step volume per minute if pump operates by speed (rpm) see page 14 'Setting the tube size'.

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Cascading several pumps

Providing that an appropriate software is available, up to 8 ISMATEC® pumps can be controlled via one PC.

In order to establish a cascade you need to connect the pumps with the connecting cable (AG0013) as shown on the left.

The same cable AG0013 serves to connect the first pump to the PC, provided that it has a 9-pin RS232 connector.

If the PC is equipped with a 25-pin connector, you need an additional adapter XC0009.

➔ **Each pump must be allocated its own individual address.**

For this procedure, you first have to connect each pump in turn to the RS232 interface of the PC. Two or more connected pumps must not have the same address. Only when all the pumps have been assigned their addresses can you cascade them as shown on the left.

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

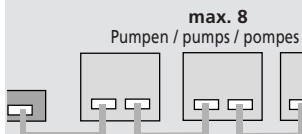
Each pump requires an individual address.
(see page 28).

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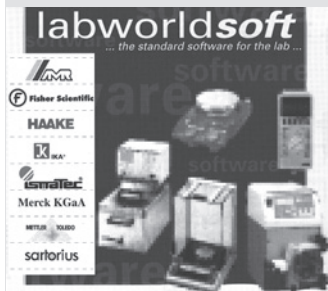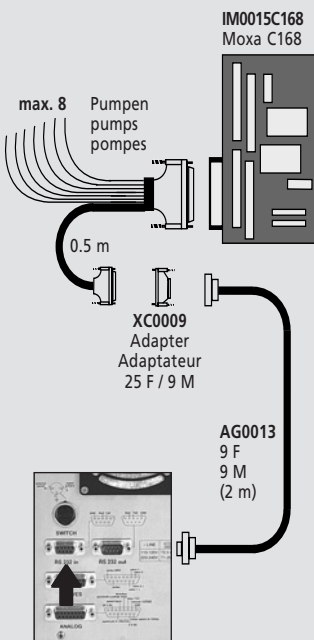Programming Software

Labworldsoft®

is an innovative Windows software for controlling, measuring and evaluating operations in connection with a number of different laboratory instruments.

Hardware requirements

486 PC, at least 33 MHz and 8 MB RAM, mouse, VGA display, monochrome, with at least 16 grey scale values

Software requirement

Windows 3.1 or higher,
Windows 95

Labworldsoft®

Order No. SOF103

Interface card Moxa C168

including a 25 pin multiple cable (male) for controlling up to 8 laboratory instruments

Order No. IM0015C168

Adaptor 25 F / 9 M

25 pins, female / 9 pins, male for IM0015C168 (1 adapter for each controlled instrument)

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Accessories

Foot switch

Order No. ISM 894
This foot switch serves as a start/stop device. It is very useful when using the pump as a dispenser for filling tubes, bottles, etc.. Both hands are free for handling the bottles and tubing.

Extension cable 9F / 9M

9 pins, female / 9 pins, male
für XC0009 (2 m long)
Order No. AG0013

Adaptor 25F / 9M

25 pins, female / 9 pins, male
Order-No. XC0009

Titration dispenser

in preparation

Spare-cassettes

Cassette MS/CA Click'n'go
Material **POM**
Order No. **IS 3510**

Cassette MS/CA pressure lever
Material **POM**
Order No. **IS 0649**

Cassette MS/CA pressure lever
Material **PVDF**
Order No. **IS 3629**

- ➔ The cassettes with pressure lever are available on request. This type of cassette may provides better results both for the Tygon MH tubing and at elevated differential pressure conditions.
- ➔ The cassettes allow the user to insert tubes with different diameters and materials on the same roller-head.

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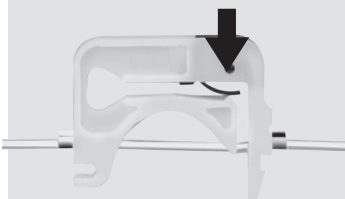
MS/CA Anpresshebel

MS/CA pressure lever
MS/CA levier de pression

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△ Please note

The values indicated in this table are only approximate and determined as follows:

ml/min, per channel, with water and Tygon® tubing, without differential pressure

Order No. = Tygon® ST (R3603)

We will be pleased to help the user to select the tubing material.

However, the user himself has the final responsibility for the selection of the correct tubing material.

For reproducible results we recommend you to use tubing from ISMATEC®.

Please ask for the detailed tubing documentation.

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Maintenance

Provided the REGLO *Digital* is operated properly and in compliance with this manual, the tubing is the only part that is subject to wear and tear.

Service and repairs

For repairs send the complete pump with detailed description of the failure to your ISMATEC® distributor.

Replacement parts

For repairs you intend to carry out by yourself (out of the warranty period), you can ask your ISMATEC® distributor for:

- parts lists
- replacement parts for pump drive
- wiring diagrams

Please give information on defect, date of purchase, serial no., and model.

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△ Please note

We also recommend you to observe our **Warranty Terms as well as our Terms and Conditions of Sale**.

In case of any queries, please contact your local ISMATEC® representative.

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2 channels 1.6–160 rpm

4 channels 1.0–100 rpm

digitally adjustable in steps of 0.1 rpm

Differential pressure

max. 1.0 bar (100 psi)

Remote control

via RS232 interface (see pages 27 to 34)

Mains connection

115/230 V_{AC} / 50/60 Hz

Fuse rating

2 x 250 mA, slow-blow (230 V_{AC})

2 x 500 mA, slow-blow (115 V_{AC})

Power consumption

max. 20 W

Protection rating

IP 30

Operating conditions

Temperature +5 to +40°C,

Rel. humidity max. 80%

- not condensing, at normal environmental conditions

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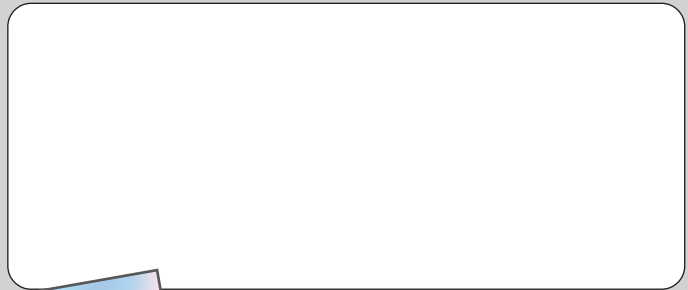
Ask for the latest
ISMATEC-
catalogue.

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ISMATEC® - Your competent partner for demanding metering and dispensing applications.

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A Unit of IDEX Corporation
Laboratoriumstechnik GmbH